

TRANSDUCER CONNECTOR INTERFACES (TCI)






- BSL-TCI0 Interface Grass
- BSL-TCI1 Interface Beckman 5 pin
- BSL-TCI2 Interface WPI, ADI, iWorx
- BSL-TCI3 Interface Lafayette
- BSL-TCI4 Interface Honeywell
- BSL-TCI5 Interface Mod Phone
- BSL-TCI6 Interface Beckman
- BSL-TCI7 Interface Nihon Koden
- BSL-TCI8 Interface Narco 7
- BSL-TCI9 Interface Fukuda
- BSL-TCI10 Interface Gould: *discontinued*
- BSL-TCI11 Interface Hugo Sachs
- BSL-TCI12 Interface Thornton

- BSL-TCI13 Interface MP30 to Piezo
- BSL-TCI14 Interface 1/4 phono
- BSL-TCI15 Interface 5-pin DIN Vernier
- BSL-TCI16 Interface BT connector, Vernier
- BSL-TCI17 Interface 5-pin Intellitool
- BSL-TCI18 Interface 2 mm Hg Strain
- BSL-TCI19 Interface 6-pin Intellitool
- BSL-TCI20 Interface 3.5 mm Intellitool
- BSL-TCI21 Interface BNC pH probe
- BSL-TCI22 Interface SS2L to SS39L

See also: SS-KIT-IN BSL/SS Custom Input Kit
SS-KIT-OUT BSL/SS Custom Output Kit

BSL-TCI PIN-OUTS

Check your hardware pin-out to match the BSL-TCI pin-out below

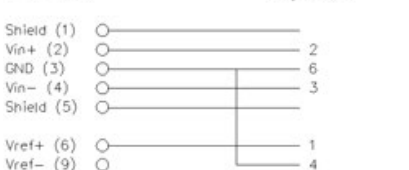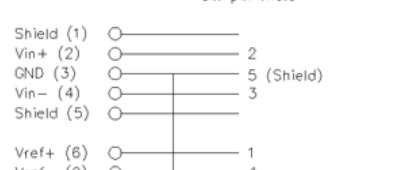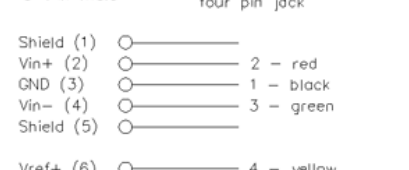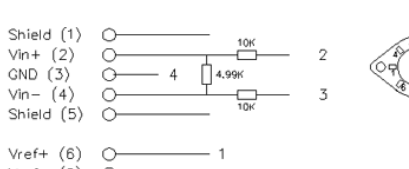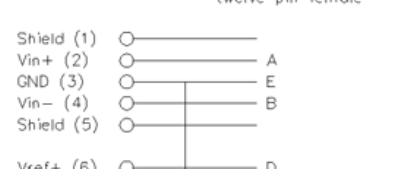BSL-TCI0 <u>BSL-TCI0 (Grass)</u> 9-Pin male Connector Pin-out, six pin male 	BSL-TCI4 <u>BSL-TCI4 (Honeywell)</u> 9-Pin male Connector Pin-out, six pin male 
BSL-TCI1 <u>BSL-TCI1 (Beckman 5-pin)</u> 9-Pin male Connector Pin-out, five pin female 	BSL-TCI5 <u>BSL-TCI5 (Modular Phone Jack)</u> 9-Pin male Connector Pin-out, four pin jack 
BSL-TCI2 <u>BSL-TCI2</u> 9-Pin male Connector Pin-out, 8-Pin female DIN 	BSL-TCI6 <u>BSL-TCI6 (Beckman 12-pin)</u> 9-Pin male Connector Pin-out, twelve pin female 


BSL-TCI PIN-OUTS

Check your hardware pin-out to match the BSL-TCI pin-out below

BSL-TCI3

BSL-TCI3 (Lafayette)


9-Pin male Connector Pin-out,
nine pin female

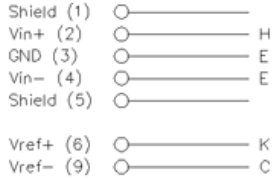

BSL-TCI7

BSL-TCI7 (Nihon Koden)


9-Pin male Connector Pin-out,
five pin female

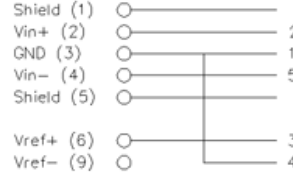

BSL-TCI8

BSL-TCI8 (Narco)


9-Pin male Connector Pin-out,
seven pin female

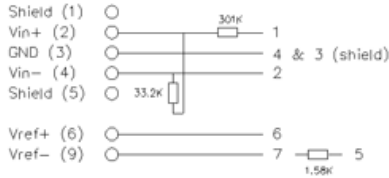

BSL-TCI14

BSL-TCI14 (Lafayette Phono)


9-Pin male Connector Pin-out,
Phono female

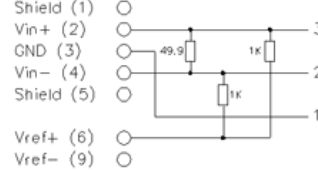

BSL-TCI9

BSL-TCI9 (Fukuda)


9-Pin male Connector Pin-out,
eight pin female

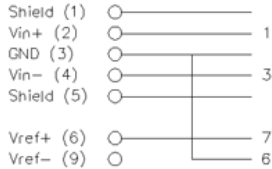

BSL-TCI15

BSL-TCI15 (Vernier 5-Pin)

9-Pin male Connector
5-PIN DIN FEMALE

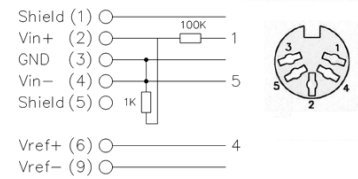


BSL-TCI10 Gould 12-pin: discontinued

BSL-TCI11

BSL-TCI11 (Hugo Sachs-Harvard)


9-Pin male Connector Pin-out,
six pin female

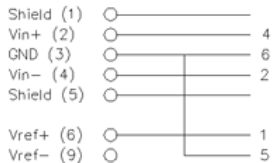

BSL-TCI16

BSL-TCI16 (Vernier Dissolved O2)


9-Pin male Connector Pin-out,
BT female

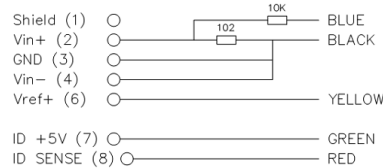

BSL-TCI12

BSL-TCI12 (Thornton)


9-Pin male Connector Pin-out,
six pin female

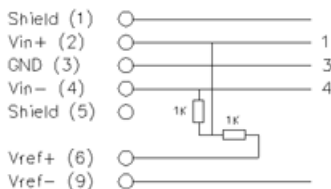

BSL-TCI17

BSL-TCI17 (Intellitool Physiogrip)

9-Pin male Connector Pin-out,
5-Pin male

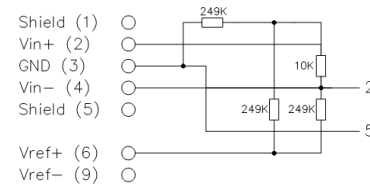


BSL-TCI PIN-OUTS

Check your hardware pin-out to match the BSL-TCI pin-out below


BSL-TCI13

BSL-TCI13 (MP3X to Piezo)

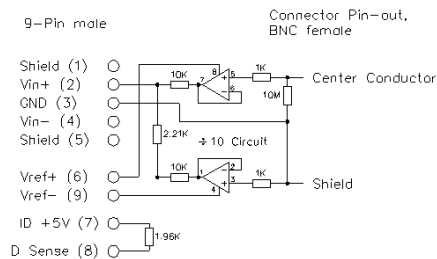


BSL-TCI18

BSL-TCI18 (Liquid Metal Strain Gauge)

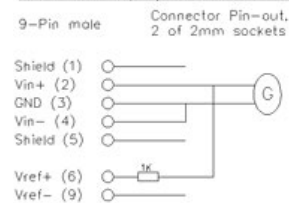


BSL-TCI19

BSL-TCI19
(Intellitool R. Hammer, DIN)

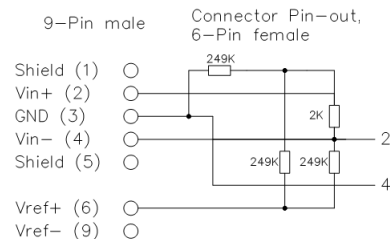


BSL-TCI21

BSL-TCI21 (BNC pH)

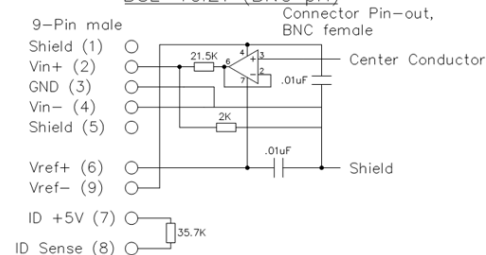

BSL-TCI20

BSL-TCI20 (Intellitool R. Hammer, phono)

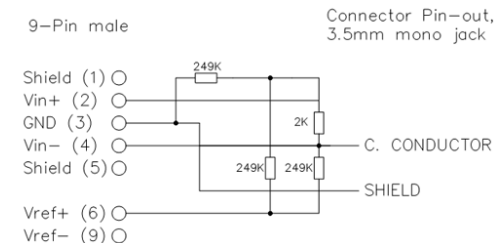

BSL-TCI22

